

USER MANUAL

T1-switch to Channel 1 timer

T2-switch to Channel 2 timer

M-set memory and return to memory timer

HR-adjust Hour

MIN-adjust Minute

SEC-adjust Second

START/STOP-start or pause the timer

HIGH/LOW/OFF slide switch-the timer alarm beep volume control

HOW TO USE

Press T1 or T2 to switch the channel the user wants to operate.

Set the timer

- 1.Press HR button to adjust “hour”, press and hold it can adjust the value faster.
- 2.Press MIN button to adjust “minute”, press and hold it can adjust the value faster.
- 3.Press SEC button to adjust “second”, press and hold it can adjust the value faster.
- 4.Press and hold HR and MIN can reset the timer to 00:00:00

The max time of each timer is 99 hour 59 min 59 sec.

Start the timer

- 1.After finishing the timer setting, press START/STOP button to start countdown timer function, the countdown function indication icon ▼ will flashes; press START/STOP again to pause it., and the icon ▼ will stop flashing; press it again, the countdown timer will continue to work.
- 2.When the countdown timer is working, the channel button green LED light indicator that the user selects will flash.

After the countdown finishes

- 1.When the timer is 00:00:00, the count-up function indication icon ▲ will flash, and the timer will start from zero. The buzzer will beep for 1 minute (the alarm beep is on), and the red LED light indicator will flash for 1 minute.
- 2.Press START/STOP once, the preset timer will show up again.

How to set the memory timer

The user may set some the same time often, the memory timer function will make this operation easier.

- 1.When the timer is not working, press and hold M button till the digits flash.
- 2.Press HR button to adjust “hour”, press and hold it can adjust the value faster.
- 3.Press MIN button to adjust “minute”, press and hold it can adjust the value faster.
- 4.Press SEC button to adjust “second”, press and hold it can adjust the value faster.
- 5.Press and hold HR and MIN can reset the timer to 00:00:00

Whenever the user wants to use the memory timer, press the M button once, and then the memory timer shows up.

How to use the stopwatch function

Press T1 or T2 to select the channel.

- 1.When the timer is paused, press and hold HR and MIN together to reset the timer.
- 2.Press START/STOP button once, the stopwatch function starts to work.

Remarks: The stopwatch function is laser tuned for 1/100th second accuracy. This stopwatch times 1/100th second for 1 hour, then 1 second increments up to 99 hour 59 min 59 sec.

Adjustable beep volume

Switch to adjust alarm beep sound (HIGH/LOW/OFF)

How to place it

Magnetic back, retractable stand and hanging bracket for hook.

How to replace the battery

When the battery power is low, the display will be dark, and need to replace new batteries with 2XAAA